


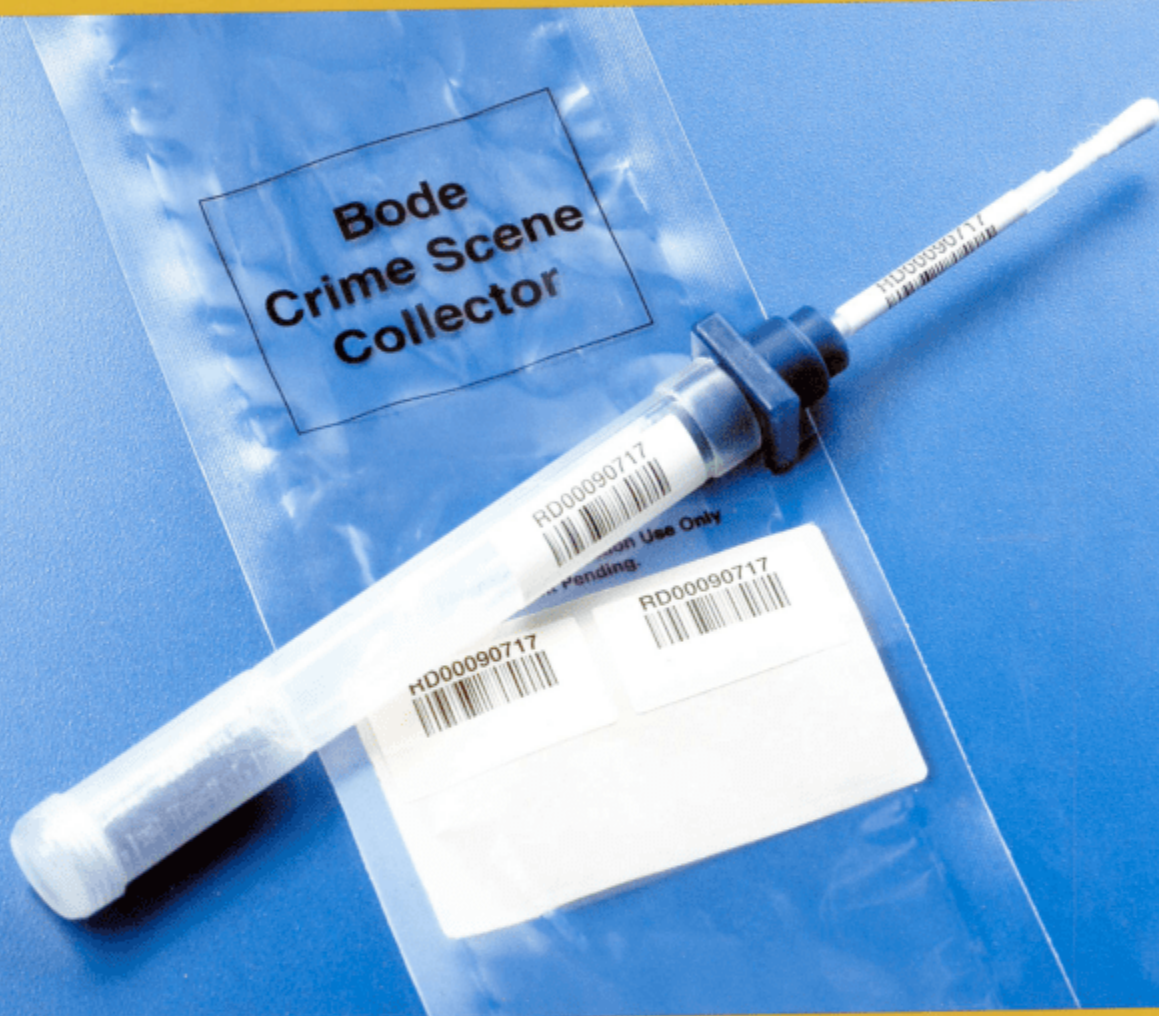

The Bode Crime Scene Collector™

The faster, more effective way to secure reliable evidence from crime scenes.

Act with confidence.

Secure evidence more efficiently, in less time, with the Bode Crime Scene Collector™



Introducing the unique Bode Crime Scene Collector™ — from the leader in DNA and advanced crime scene technologies and services.

The Bode Crime Scene Collector makes evidence collection faster and easier.

You just swab, seal, and ship. The collector maintains chain of custody, ensures evidence stability and provides protection of evidence. Components are pre-barcoded and include matching barcode labels for chain of custody of documents. The innovative shaft design prevents the shaft from breaking during the collection process and simplifies the removal of the cotton swab for processing. Evidence can be air dried or dried in the protective tube with the indicating desiccant. The collector can also be configured into kits with easy-to-follow instructions and all materials required for simple collection.

The indicating desiccant is designed to dry a sample in 16 hours when used as directed. A box of 50 collectors is packaged in five foil pouches, each containing 10 collectors and 10 Crime Scene Transport Pouches. The foil pouch protects the indicating desiccant from moisture until opened. The indicating desiccant will change color upon saturation and can be replaced if needed.

Features

- Package protects collector from contamination and moisture before use.
- Integrated indicating desiccant dries evidence to ensure stability during transportation to crime lab.
- Collectors are pre-labeled with unique barcodes — no two collectors have the same number.
- Additional matching barcodes are on collector package for photo documentation of evidence or labeling crime scene documents and transport packages.
- Traditional cotton-tipped swab allows for wet or dry evidence collection.
- Ergonomic plastic tube/handle easily converts into protective, secure, polypropylene transportation device in one motion.
- Collector design minimizes contamination of evidence, errors in chain of custody and processing time.

A complete evidence collection system in one easy tool.



- A. cotton swab**
- B. reversible cap**
- C. protective tube with barcode label**
- D. indicating desiccant**


Order today by calling 866.BODE.4.ID extension 787.

Product Catalog Numbers: Bode Crime Scene Collector P01D58 (box of 50) and P01D59 (case of 4 boxes).




COLLECTING

The Bode Crime Scene Collector is faster and more reliable than traditional methods because it's easier to open, hold, use and reseal. The pre-labeled collector and additional barcodes also reduce labeling time.

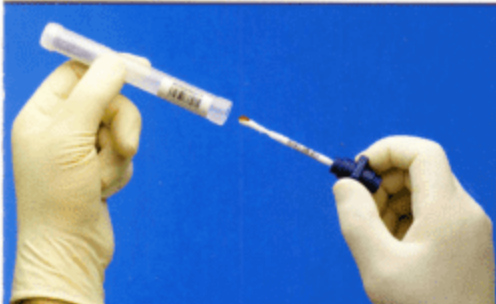

PROTECTING

Inserting the swab into the protective tube and placing the collector into the Crime Scene Transport Pouch secures evidence in a sealed environment.

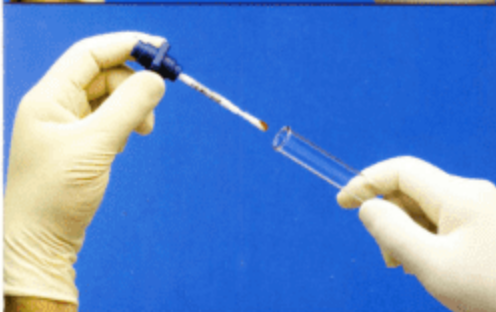

PROCESSING

The swab is easily transferred from the collector for automated processing, minimizing errors and contamination. Just hold the Crime Scene Collector in a tube at a 45-degree angle and apply pressure to snap off the cotton swab.

INSTRUCTIONS

Step 1

Place collector wrapper with visible barcode number near evidence for photo documentation. Remove wrapper barcode labels for identifying crime scene documents and labeling outer package.

Step 2

Tear open wrapper and remove collector.

Step 3

For dry stain collection, add 100 microliters of sterile water (about 3 drops) to the swab then touch and rotate the swab on stain. For wet evidence, dab the swab in the evidence to collect about 100 microliters. Air drying will be necessary if too much evidence is collected.

Step 4

Firmly grip cap and pull protective tube away.

Step 5

Turn cap 180° and insert swab into tube. Seal by pressing tube firmly onto cap.

Step 6

Insert collector into paper transport pouch and seal. Remove a barcode label and apply to transport pouch for identification.

SPECIFICATIONS

Length

8 3/8", converts to 6" transportation tube after evidence collection

Barcodes

128C symbology with two alpha and eight numeric digits

Desiccant

Two-gram capsule in bottom of tube/handle; desiccant turns color as it's saturated

Drying Time

Three drops of liquid on swab dries overnight

Packaging

50 collectors and 50 Crime Scene Transport Pouches per box, packaged in five resealable foil pouches